

IGOR template: Import tools for MED Conductor data

1. ReadData.ipf: Data import from all 64 channel & one sweep.
(Command) OpenMEDFile(nTargetSw)
(Argument) nTargetSw: sweep number to focus
(Other operation) Select the data file name in the file dialog window during the program.
(Result) This command will load the data of each channel into wave format, "ch#" (# is number of channel).

2. ReadDataOneCh.ipf: Data import from one channel & all sweep.
(Command) OpenMEDAllSw1ch(nTargetCh)
(Argument) nTargetCh: channel number to focus
(Other operation) Select the data file name in the file dialog window during the program.
(Result) This command will load the data of each sweep into wave format, "sw#" (# is number of sweep).