

MED Wrapper Excel Macro


1. Installation of the DLL file in your PC

In order to use the function in this wrapper, you have to installation in the Windows and Excel.

- 1) Copy “MedWrapper.dll” in the somewhere in PC.
(e.g. C:\Medwrapper).
- 2) Run MS-DOS prompt and move to the holder, where you copy the dll file
(e.g. > CD C:\MedWrapper)
- 3) Install to the registry in the windows.

In the case of Windows 2000: > C:\windows\system32\regsvr32.exe MedWrapper.dll

If installation will succeed, the following window pops up.

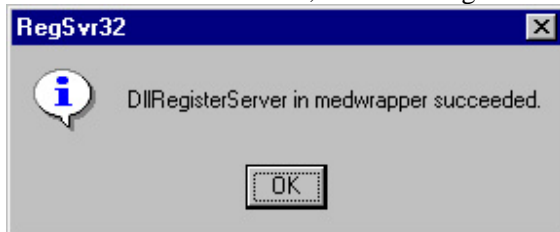

- 4) Install to the Excel
 - a) Run the “MED_Conductor_Macro.xls”, and change “Visual Basic Window” with pushing [Alt]+[F11]
 - b) Select the [tool] – [references] in the menu bar.
 (“References- VBA project” window should pop up.)
 - c) Push [Browse...] button and select the “MedWrapper.dll”, which you saved before.
 - d) Save it
 - e) Close the window

NOTE: if you would doubt this setting, please check the following procedures.

- a) In “Visual Basic Window”, select [view] – [object browser] in the menu bar.
- b) “Object browser” window would pop up
- c) If you select “MedWrapper” at “<All Libraries>” window on the left upper side, “Meddata” should appear in the class window. If not, could you check again above procedures.